

INSTRUCTION FOR USE

P.A.S. Periodic Acid Schiff

<u>Catalog number</u>	<u>Size</u>
04-130802	100 test

Expected aim

Product for the preparation of cyto-histological samples for optical microscopy.

Method :

- 1) Method for histological sections
 - Bring section to distilled water.
 - Put on the section 10 drops of reagent A : leave to act 10 minutes.
 - Wash in distilled water.
 - Put on the section 10 drops of reagent B : leave to act 20 minutes.
 - Wash in distilled water.
 - Put on the section 10 drops of reagent C: leave to act 2 minutes.
 - Drain the slide and without washing add 10 drops of reagent D: leave to act 2 minutes. Rinse in distilled water.
 - Put on the section 10 drops of reagent E: 3 minutes.
 - Wash in running tap water for 5 minutes.
 - Dehydrate through ascending alcohols. Clear in xylene and mount.

- 2) Method for hematology and cytology
 - Bring section to distilled water.
 - Put on the smear 10 drops of reagent A : leave to act 10 minutes.
 - Wash in distilled water.
 - Put on the smear 10 drops of reagent B : leave to act 15 minutes.
 - Wash in distilled water.
 - Put on the smear 10 drops of reagent C: leave to act 2 minutes.
 - Drain the slide and without washing add 10 drops of reagent D: leave to act 2 minutes. Rinse in distilled water
 - Put on the smear 10 drops of reagent E: 3 minutes.
 - Wash in running tap water for 5 minutes.
 - Dehydrate through ascending alcohols. Clear in xylene mount.

Warning and Precaution

The product must be used exclusively by specialized technical operators.
Carefully read the information on the classification of dangerous substances on the label. Always refer to the safety data sheet where are available the information on the risk presented by the mixture, the precautionary measures during use, the measures first aid and the event of accidental release.
Do not use if the primary container is damaged.

Contraindication

Obtain special instructions before use.
Wear protective gloves / clothing and eye / face protection.
Keep away from heat/sparks/open flames/hot surface. No Smoking.
IF ON SKIN (on hair): remove/take off immediately all contaminated clothing. Rinse skin with water/shower.