

INSTRUCTION FOR USE

Silver Impregnation

<u>Catalog number</u>	<u>Size</u>
04-040801	100 test

Expected aim

Product for the preparation of cyto-histological samples for optical microscopy.

Method

- 1) Bring the section to distilled water.
- 2) Put on the section 5 drops of reagent A and 5 drops of reagent B: leave to act 5 minutes.
- 3) Wash the slide in distilled water.
- 4) Put on the section 10 drops of reagent C: leave to act 3 minutes.
- 5) Double washing in distilled water.
- 6) Put on the section 10 drops of reagent D: leave to act 3 minutes.
- 7) Double washing in distilled water.
- 8) Put on the section 10 drops of reagent E: leave to act 3 minutes.
- 9) Wash in distilled water.
- 10) Put on the section 10 drops of reagent F: leave to act 5 minutes.
- 11) Double washing in distilled water.
- 12) Put on the section 10 drops of reagent G: leave to act 5 minutes.
- 13) Wash in tap water 5 minutes.
- 14) Dehydrate through ascending alcohols: clear in xylene and mount.

Warning and Precaution

The product must be used exclusively by specialized technical operators.

Carefully read the information on the classification of dangerous substances on the label. Always refer to the safety data sheet where are available the information on the risk presented by the mixture, the precautionary measures during use, the measures first aid and the event of accidental release. Do not use if the primary container is damaged.

Contraindication

Obtain special instructions before use.

Wear protective gloves / clothing and eye / face protection.

Keep away from heat/sparks/open flames/hot surface. No Smoking.

IF ON SKIN (on hair): remove/take off immediately all contaminated clothing. Rinse skin with water/shower.